

ATCC Medium 1445: Nutrient Agar/Broth w/ Streptomycin 20µg/mL

Agar Medium:

Nutrient Agar..... 23.0 g
DI Water..... 950.0 mL

Broth Medium:

Nutrient Broth..... 8.0 g
DI Water..... 950.0 mL

Mix media well and autoclave at 121°C.
Cool to 45-50°C and add the following solution:

Streptomycin Solution

Streptomycin..... 20.0 mg
DI Water..... 50.0 mL

Filter Sterilize.